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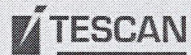
Conditions:

Sample Gold-on-carbon AGAR S1987 at position 7 (center) of the sample carousel.

Specimen AGS142 for STEM detector.

Chamber and resolution test sample decontaminated. Recommended time is: 30min without any sample in the chamber and then 20s (not longer, stop it manually if it is running longer than 20s) with resolution test sample with stage in Home position.

Description of the test:

1. All images should be saved with following format: 1024 x 768 pixels, 16bit grayscale, PNG format.
2. Scanning speed: 6.
3. Line accumulation: 4x.
4. BC optimal (Set Best Resolution) or as specified by the table below.
5. Magnification: 300kx and 1000kx. Images at 300kx are used for assessing general image quality by eye. Images at 1000kx are used for resolution measurement.
6. For BDM mode imaging use always position 7 of the sample carousel.

Detector	Mode	WD [mm]	HV [kV]	Beam Current	Magnifica tion	Image name	Typical Resolution [nm]	Measured Resolution [nm]
In-Beam SE	UH- Resolution CrossFree mode	3.5	15	CrossFree	1000kx	RT_IBSE_15_1000	0.6+10%	
					300kx	RT_IBSE_15_300	N/A	N/A
In-Beam SE	UH- Resolution	1	1	5 pA	1000kx	RT_IBSE_01_1000	1.2+10%	
					300kx	RT_IBSE_01_300	N/A	N/A
In-Beam BSE	UH Resolution	3.5	15	100 pA	1000kx	RT_IBBSE_15_1000	1.4+10%	
					300kx	RT_IBBSE_15_300	N/A	N/A
SE (BDM mode)	UH Resolution CrossFree mode	2	1	CrossFree	1000kx	RT_BDM_05_1000	0,9+10%	
					300kx	RT_BDM_05_300	N/A	N/A

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1. Turn ON the FIB emission (2 μ A).
2. After emission is running, wait for 5 minutes.
3. After 5 minutes note the value of Suppressor voltage (Initial value).
4. Wait for 1 hour and note the Suppressor voltage value (After 1 h value). If emission is stable, the Suppressor voltage should not change more than 500 V.

Suppressor initial value [kV]	
Suppressor value after 1 h [kV]	
Difference [V]	

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Conditions:

- Use silicon wafer at FIB-SEM intersection.
- FIB emission stable and repeatable.
- SEM observation parameters: 10 kV, 300 pA, DEPTH

Description of the test:

1. Check all factory presets.
2. If necessary, adjust FIB probe current to be within 5 % of specified current of the preset by changing the value of FIB Condenser voltage. The preset 1 pA must be ≤ 1 pA and the preset 100 nA must be ≥ 100 nA.
3. Make sure that there is no image movement on FIB OBJ wobble for presets less than 500 pA.
4. Measure the milled spots according to the table.
5. Save re-tuned presets, if necessary.
6. Save acquired images of spots into Acceptance folder.

ORAGE at HV 30 kV – milling for 3 sec		
FIB beam current	Specified hole diameter	Measured hole diameter
100 nA	$\leq 11 \mu\text{m}$	
50 nA	$\leq 6.5 \mu\text{m}$	
1 nA	$\leq 0.9 \mu\text{m}$	
50 pA	$\leq 0.3 \mu\text{m}$	

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Conditions:

- FIB Preset: 30 keV; UHR imaging; E-T detector.
- Sample Agar S1937 (tin-on-carbon).
- Sample tilt 55° ; sample in the intersection point.
- FIB perfectly centered (including aperture).
- Image parameters: resolution 1024 x 1024 pixels, bit depth: 16 bits, format PNG.
- FIB scan speed: 6.
- Field of view $1 \mu\text{m}$.

Description of the test:

1. Acquire image and save it as: RT_FIB.png.

- Use ResolutionGUI tool for R50 resolution measurement. Measured value must be less or equal to the specified value.

FIB resolution	Measured [nm]
Typical value: $\leq 2.5_{+10\%}$ [nm]	

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Conditions:

- Use plain Si wafer for tests.
- Heating of tested precursors should be on for at least 15 minutes.
- Sample at SEM-FIB intersection, sample tilt = 55 deg
- FIB acceleration voltage: 30 keV.
- FIB preset 30 keV, 250 pA
- View field: $\approx 25 \mu\text{m}$

Description of the test:

- Run Draw Beam, project Protection_layer_deposition, do not change the project settings it should be correct (Spot size = 50nm, Spacing = 2, Dwell time = 0.1 μs , Scanning Path: Zig-zag, Exposure pitch = 100 nm, Exposure time = 180 s, rectangle size 10x1 μm)
- Choose the precursor and start the deposition. Check the chamber pressure stability during the deposition. The pressure should be stable and change rate less than $1 \cdot 10^{-4}$ Pa/3min. If the condition is not met, repeat the deposition
- Choose: SEM \rightarrow Geometric transformation \rightarrow Tilt corr. mode: follow sample cross section (FCS) for deposition of thickness measurement.
- Measure the height and save the image of the deposited rectangle.

Precursor	FIB beam current used [pA]	Minimal deposited height [μm]	Measured deposition height [μm]
Platinum		>1	
Carbon		>1	

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Conditions:

- All GIS in parking positions.
- Room temperature stability better than 2°C with the rate of $1^\circ\text{C}/\text{hour}$ ($0.017^\circ\text{C}/\text{min}$).
- Calibrate all nanomanipulators before test.
- SEM Imaging conditions: 5 kV.
- FIB Imaging conditions: 30 keV HR Imaging.
- No sample in SEM-FIB intersection.

Description of the test – repeat the test for each nanomanipulator arm:

1. Move nanomanipulator arm into intersection, you should see nanomanipulator needle in both SEM and FIB scanning windows. Save the position into the memory.
2. Change the scanning speed to 5 and re-scan both FIB and SEM images. Mark position of the manipulator needle in both FIB and SEM window.
3. Move nanomanipulator to parking position and back to the saved intersection position.
4. Re-Scan both FIB and SEM image, and measure nanomanipulator needle shifts in both FIB and SEM. Write down the difference into Work-park-work row. The maximum allowed difference is 5 μm .
5. Mark position of the needle again and wait for 5 minutes, write down the needle drift within 5 minutes into the table. Maximum allowed drift is 5 μm

	SEM scan window	FIB scan window
Work-park-work [μm]		
5 min needle drift [μm]		
Max. differences [μm]	$\leq 5 \mu\text{m}$	$\leq 5 \mu\text{m}$

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